


Environmental Allergen Swab Collection & Shipping Procedure

Sampling/Swabbing

1. Pre-moisten sterile swab by dipping the swab in the extraction solution provided.
2. Using a sterile swab, sample an area 4 in. x 4 in. using a crosshatch technique.
3. Place swab into original tube once sampling is complete. Remember to label each tube.

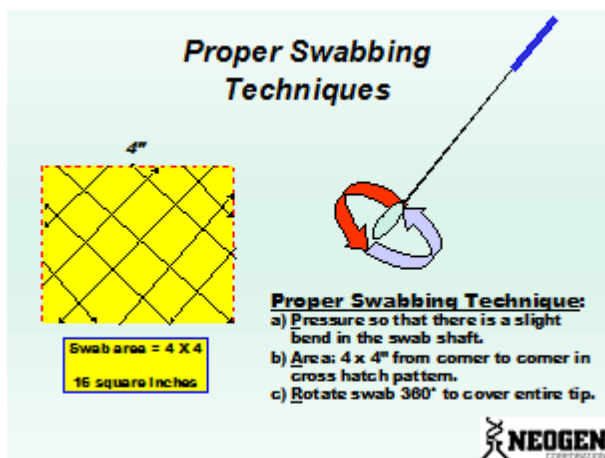

Preparing Swab for Shipping

Preparing Swab for Shipping

Add a few drops of extraction solution to the bottom of the tube (do not fill higher than the bottom of the swab bud—see arrow at left).

Replace swab in tube, making sure it is sealed tightly. Store swabs in the refrigerator (2–8°C) until ready to ship.

Ship prepared swabs to their final destination using overnight service. Include an ice pack with the swabs to keep them cool during shipping.

NOTE: Swabs should arrive at the testing destination within 24 hours of collection.

*Note please do not submerge the swab as this invalidates the test

